

**Product Code . EL-A-11063**

## Table Top Front Loading Autoclave



### Description

---

#### Table Top Front Loading Autoclave

##### Features:

Auto drain of water to reservoir tank after cycle completion

Completely automatic & with error message

Table Top Front loading Sterilizer / Autoclave

Digital Pressure Display & also Analogue pressure display

Seamless chamber

Feasibility to take out instruments from chamber even after electricity fails

Fast cycle time with Digital timer for Wet & Dry Cycle

##### Mode Select:

Liquid programs to sterilize liquid sterilization temperature : 121°C, sterilization time : 20 minutes & dry time : 15 minutes.

Instrument programs to sterilize metallic devices sterilization temperature : 134°C, sterilization time :

---

4 minutes & dry time : 15 minutes.




Rubber programs to sterilize rubber & plastic sterilization temperature : 121°C, sterilization time : 20 minutes & dry time : 15 minutes.

### Manual Program:

Manual Program Manual program as per user requirements program can set pre vacuum, post vacuum sterilization temperature : 121°C~134°C, sterilization time : 1~60 minutes, dry time : 1~60 minutes.

```
{ "@context": "https://schema.org/", "@type": "Product", "name": "Table Top Front Loading Autoclave", "image": "http://www.educational-equipments.com/images/catalog/product/848083597TableTopFrontLoadingAutoclave.jpg", "description": "Table Top Front Loading Autoclave Features: Auto drain of water to reservoir tank after cycle completion Completely automatic & with error message Table Top Front loading Sterilizer / Autoclave Digital Pressure Display & also Analogue pressure display Seamless chamber Feasibility to take out instruments from chamber even after electricity fails Fast cycle time with Digital timer for Wet & Dry Cycle Mode Select: Liquid programs to sterilize liquid sterilization temperature : 121°C, sterilization time : 20 minutes & dry time : 15 minutes. Instrument programs to sterilize metallic devices sterilization temperature : 134°C, sterilization time : 4 minutes & dry time : 15 minutes. Rubber programs to sterilize rubber & plastic sterilization temperature : 121°C, sterilization time : 20 minutes & dry time : 15 minutes.", "brand": "Educational Lab Equipments", "sku": "5", "gtin8": "5", "gtin13": "5", "gtin14": "5", "mpn": "5", "aggregateRating": { "@type": "AggregateRating", "ratingValue": "5", "bestRating": "5", "worstRating": "0", "ratingCount": "10" } }
```

---

Educational Lab Equipments,  
#449, HSIIDC, Industrial Area, Saha, Haryana  
Direct Contact Details  +91-98173-19615  sales@educational-equipments.com  
 www.educational-equipments.com